

SWING

beamline

Biophysics

ORGANISMS



TISSUES



CELLS



ORGANELLES



COMPLEXES



PROTEINS



ATOMS





SAMPLES

TYPES

- Liquids, liquid crystals, lipids, (rheo)gels
- Colloids, micelles, liposomes
- Heterogeneous solutions
- Biological tissues (fibres, bones, tendons, muscles, hair)
- Bio-inspired materials

ENVIRONMENTS

- Multi Capillaries Holder (-20 to 120 °C)
- Circulation capillary holder (10 to 60°C)
- Liquid autosampler (duty cycle: 4 min)
- Multiple Gels/Solids Holder (10 to 60°C)
- Biologic Stop Flow mixer SFM400
- Traction Cell (1 to 100 N, res: 0.05 N)
- Fluigent Microfluidic MFCS-EZ system
- Linkam Temperature Stage (-196 to 600 °C)
- Anton Paar Rheometer MCR 501
- Your own environment !

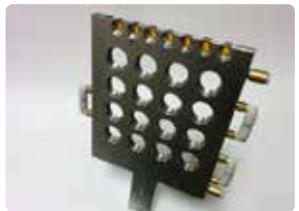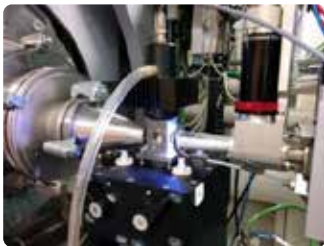

TECHNIQUE

SWING is a beamline for Small-Angle X-ray Scattering, providing structural information from supramolecular assemblies (size, shape, distances, orientations), at a scale between nm and μm . Widely used in the field of Soft Condensed Matter, it usually needs a priori models for data interpretation.



SAXS



Q-range : from $5 \cdot 10^{-4} \text{ \AA}^{-1}$ @ 6 keV
to 3.0 \AA^{-1} @ 16 keV



Typical times for one image exposure :
10 – 1000 ms.



Fast measurements down to 1.5 ms rate,
in shutterless mode.



Downloadable graphical application
for data reduction.



Typical size of one SAXS image:
1 M / 4 M pixels

μ -SAXS Cartography



Beam size (lateral spatial resolution):
10 (V) x 20 (H) μm




Continuous raster scan mode with
no dead time.



Cartography data split into interconnected
HDF5 2Gb files

HIGHLIGHTS


Designing food-grade hydrocolloidal encapsulation systems.

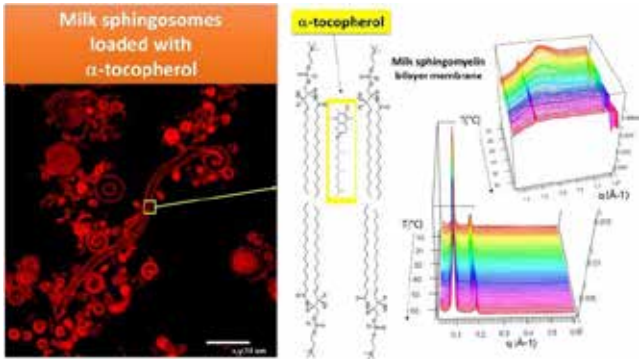

Lopez, Christelle, *et al.*, Food Research International (2022)

- Milk-sphingomyelin sphingosomes loaded with tocopherols prevent oxidation in aqueous foods containing polyunsaturated lipids such as oil-in-water emulsions.


Expanding plant-based bio-composites for structural applications, especially using flax due to its high potential.

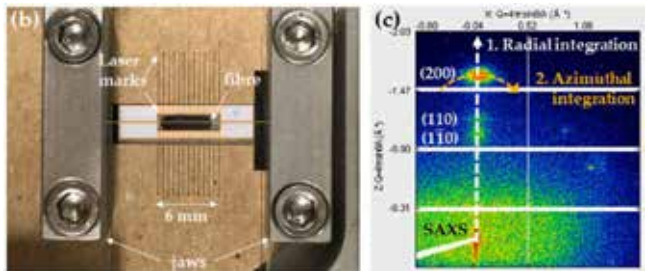

Emmanuelle Richely, Alain Bourmaud, Johnny Beaugrand *et al.*, Composites Part C: Open Access 9 (2022)

- Contrasted initial microfibril angles between 4.7 and 7.4° depending on the fibre.
- Major influence of the cellulose microfibrils on the tensile response, with a non-linear decrease of the overall Young's modulus upon increasing microfibril angle.

Hydrogels are able to retain a large amount of water content, which is of utmost importance for medicine and more generally for hygiene and medical product development.

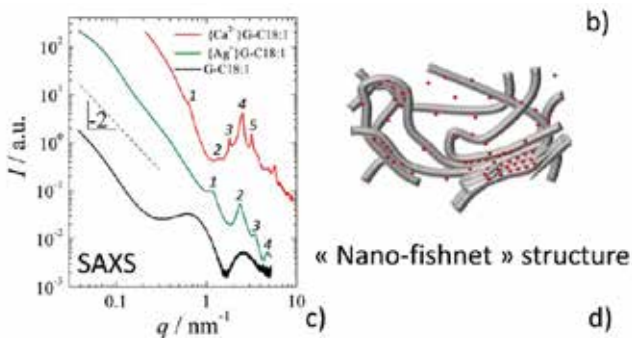

Alexandre Poirier, Niki Baccile, *et al.*, *Soft Matter* (2023)

- Strong stability of the hydrogel towards shear and temperature explained by its β -sheet-like structure.
- Unique self-assembly behaviour in the presence of Ca^{2+} or Ag^{+} , as phases from molecules with similar chemical structures are not affected by cations.

REFERENCES

- Thureau, A., Roblin, P., Perez, J. "**BioSAXS on the SWING beamline at Synchrotron SOLEIL**" *Journal of Applied Crystallography*. (2021).
- Desjardins, K., Pomorski, M., Bizien, T., Thureau, A., Meneglier, C., Pérez, J. "**An active x-ray beamstop based on single crystal CVD diamond at beamline SWING**" *Review of Scientific Instruments*. (2021).



**More information
on SWING web page**

CONTACT

Javier Pérez

Physicist, Beamline Manager

✉ javier.perez@synchrotron-soleil.fr

☎ +33 (0)1 69 35 96 19

Aurélien Thureau

Biologist, Beamline scientist

✉ aurelien.thureau@synchrotron-soleil.fr

☎ +33 (0)1 69 35 97 83

Thomas Bizien

Chemist, Beamline scientist

✉ thomas.bizien@synchrotron-soleil.fr

☎ +33 (0)1 69 35 96 20

Health & Well-Being at SOLEIL



Link to the
web page

SOLEIL's Health and Well-being Scientific Section is composed of 30 scientific experts from different fields. Through collaborative and science-driven approaches, the Section offers the community a coherent portfolio of state-of-the-art techniques to serve scientific and societal health-related challenges.



L'Orme des Merisiers - Départementale 128
91190 Saint-Aubin - FRANCE
www.synchrotron-soleil.fr